

PIPETTOR


Easy Tray Adaptor

Part Number 54563

Tray Overview

Populate 384 well plates with a 96 channel pipettor using this “Easy Tray” adaptor. It is accurate, intuitive and easy to use, a knob rotates four times to align all wells of a 384 plate under the 96 channel pipette tips. Each position positively clicks in place without needing to align by hand; the microplate is not handled during the process, eliminating the possibility of contamination.


Place a 384 well microplate into the Easy Tray. Move the knob to **Position #1** and dispense.

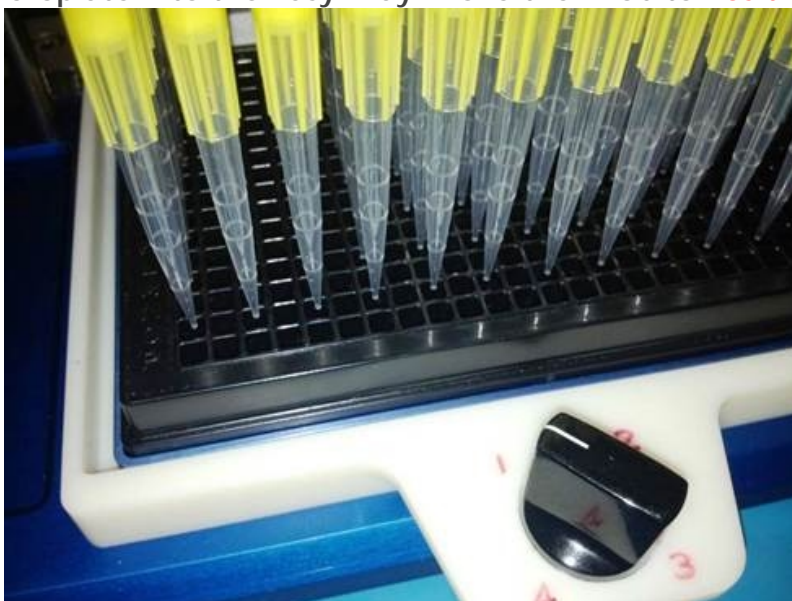


Now rotate the knob to **Position #2**, the microplate will then move to the adjacent empty well below the 96 channel tips to dispense into. Each position positively clicks in place without needing to align by hand.



Again rotate the knob to **Position #3** and the plate is positioned in the well below. Each position positively clicks in place.

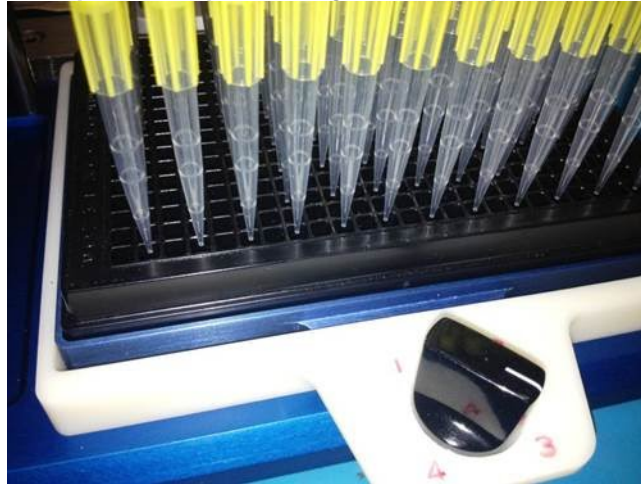

Finally rotate to **Position #4** to finish the 96 to 384 dispenses.

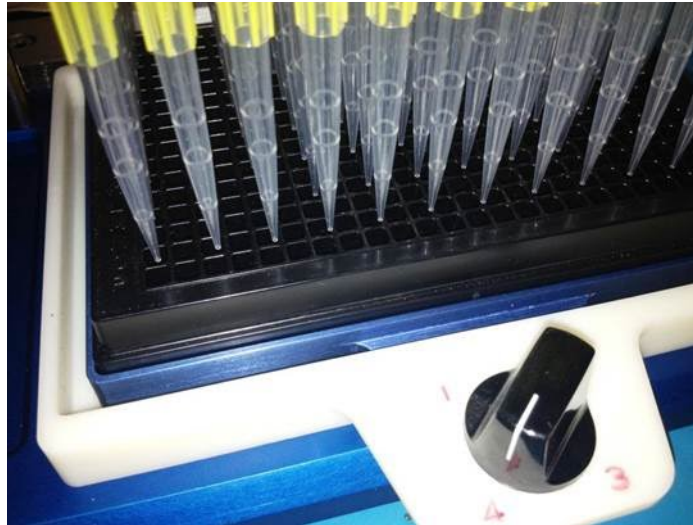

ORDER INFORMATION

Contact your local distributor
or email info@sequencebiotech.com
or contact Sequence BioTech, LLC Dan Seguin
614 Nashua Street, Unit 223
Milford NH 03055-4917
M: (603)769-0963
F: (603)732-0327